

Setting the PRBFS10-ST/PRBFS10-ST82 Probe Height for App 2

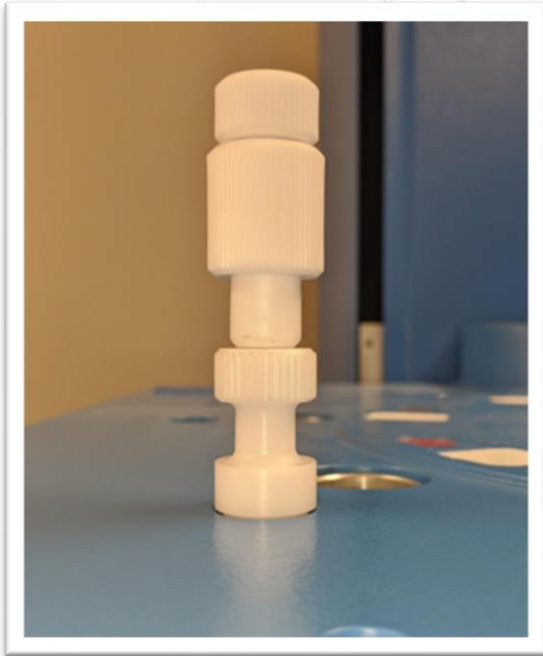


Figure 1

500 mL Sampling Position

1. Insert the PRBFS10-ST/PRBFS10-ST82 probe assembly into the position to be tested.
2. Slide the top of the probe assembly down so it is touching the bottom of the probe assembly. See Figure 1.

900 mL Sampling Position

1. Insert the PRBFS10-ST/PRBFS10-ST82 probe assembly into the position to be tested.
2. Slide the top of the probe assembly up and then insert the spacer 010-06-202.
3. Ensure the spacer is touching the top and bottom probe assemblies. See Figure 2.
4. Remove the spacer 010-06-202.



Figure 2